



System Requirements - Valab® Server V15.xx (excluding 15.01)



Characteristic	Possible values
Usage	Operation 24/7 (dedicated machine recommended, hibernation and network card sleep mode disabled). It communicates with your LIS through a TCP / IP network connection. Your LIS must have an option allowing Valab® connection and integration.
OS ⁽¹⁾	<ul style="list-style-type: none"> • Windows Server 2022® x64 Standard Edition with Desktop Experience (graphical interface) • Windows Server 2019® x64 Standard Edition with Desktop Experience (graphical interface) • Windows Server 2016® x64 Standard Edition with Desktop Experience (graphical interface) • Windows 11® Pro Edition • Windows 10® x86 and x64 Pro Edition
Processor	4-core - 64 bits - 2 Ghz
RAM	4 Gbytes
Disk type	Mechanical hard disk drive > 10 000 rpm or SSD
Typical recommended disk space	Space needed for the OS + : - licence 1200 reports / day = 50 Gbytes - licence 2400 reports / day = 75 Gbytes - licence 4000 reports / day = 110 Gbytes - licence 6500 reports / day = 170 Gbytes - licence 10 000 reports / day = 250 Gbytes - licence 15 000 reports / day = 440 Gbytes
.NET Framework	Minimum 4.5.1 version.
DB	No database engine installed
FireWall - Network	<ul style="list-style-type: none"> • Fixed IP mandatory • Authorise incoming TCP protocol connections on the port(s) used to communicate with the LIS (by default port 7200 if only one LIS connection) • Authorise incoming TCP and UDP protocol connections on the ports used to communicate with the Valab® services of the Valab® Server and with the Valab® clients (by default 8080, 8089, 8090 and 2110) • Authorise outgoing HTTPS protocol connections on the port 443 to communicate with the VSP (remote licence manager)
Access	<ul style="list-style-type: none"> • Local administrator account dedicated to the VALAB company • Remote maintenance access (VPN + RDP, remote control software,...) • Internet access to the "valab.com" domain
Backup / Security	<ul style="list-style-type: none"> • Policy and implementation to be determined by the IS Manager of the medical laboratory • We recommend that you encrypt the connection between your LIS and Valab, in particular in the case of remote hosting of these systems.



System Requirements - Valab® Client V15.xx (excluding 15.01)

Characteristic	Possible values
Usage	Thick client for remote access to the Valab® server
OS ⁽¹⁾	<ul style="list-style-type: none"> • Windows 11® • Windows 10® • Windows Server 2022® x64 • Windows Server 2019® x64 • Windows Server 2016® x64
Processor	Dual core - 64 bits - 2 Ghz
RAM	4 Gbytes
Disk space	Space needed for the OS + 20 Gbytes
Graphics card	1366 * 768 minimum recommended
.NET Framework	Minimum 4.5.1 version
FireWall - Network	Open ports 2110, 8089, 8090 and 8080 to the Valab® server
Access	Read / write access to the installation directory of the Valab® client for the user

⁽¹⁾ Subject to effective application of the latest security updates provided by Microsoft® for the OS
On condition that the OS is still maintained by Microsoft®